

# SAFETY FEATURES

- x To prevent entrapment the buddy -system must be used.
- x Two individuals must be present each time the bulk sterilizer is started.
- x The inside of the autoclave is observed to confirm that no one is inside prior to activating .
- x Both individuals will remain outside the autoclave to ensure that no one is inside as the door is closed and the sterilizer cycle is being activated.
- x By design the door to the sterilizer only closes very slowly to reduce the possibility of entrapment.
- x Emergency Shut -Off (red button) located on both sides of the autoclave.
- x